


Document Name		VisualSpreadsheet 5 How to install software	
File Name		TCS1158B VisualSpreadsheet 5 How to install software	
Code	TCS1158	Date	3/19/21
Author	K. Johnson	Reviewed By	S. Barton
Locations	K:\Customer Support\2 - Support\VisualSpreadsheet\5.X		
<b>Revision History</b>			
<b>Issue Date</b>	<b>Rev</b>	<b>Author</b>	<b>Changes</b>
7/10/19	A	KC, HAW	initial release
4/7/2020	B	KRJ	review and converted into controlled version
3/19/21	C	KRJ	added note at the end of the document re:anti-virus software concern


**Purpose:** How to install VisualSpreadsheet 5

There are a number of steps in order to install VisualSpreadsheet 5:


1. Click on the link provided to you by Fluid Imaging. The software will prepare to install onto your system. You will then see the following welcome screen. **Click NEXT.**

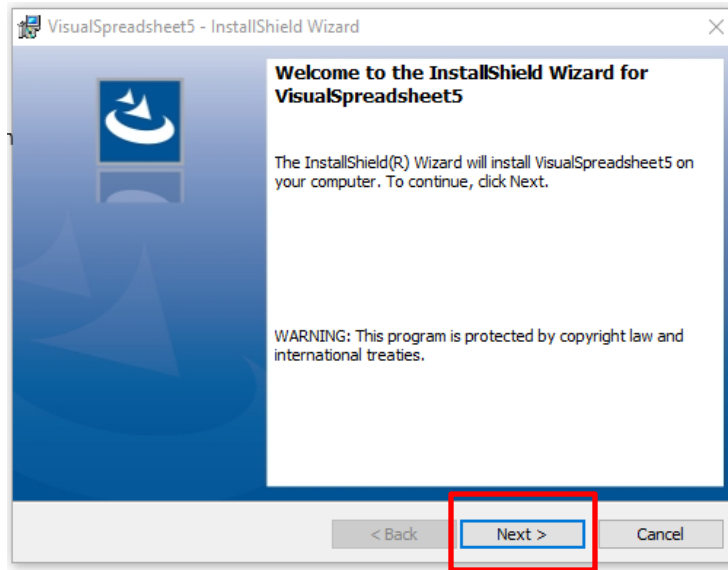


2. You will then see the End-User License Agreement. Once you have read it, please select "I accept the terms in the license agreement."

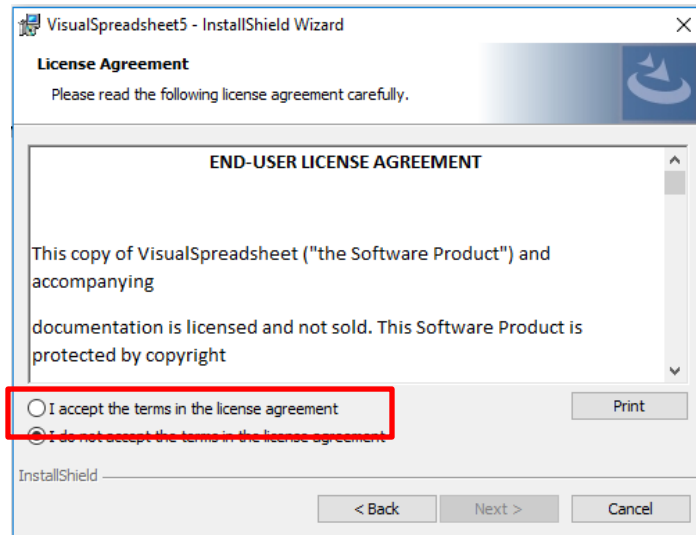


3. You will then see the Destination Folder dialog, **click NEXT**. *DO NOT* change the location of the folder.

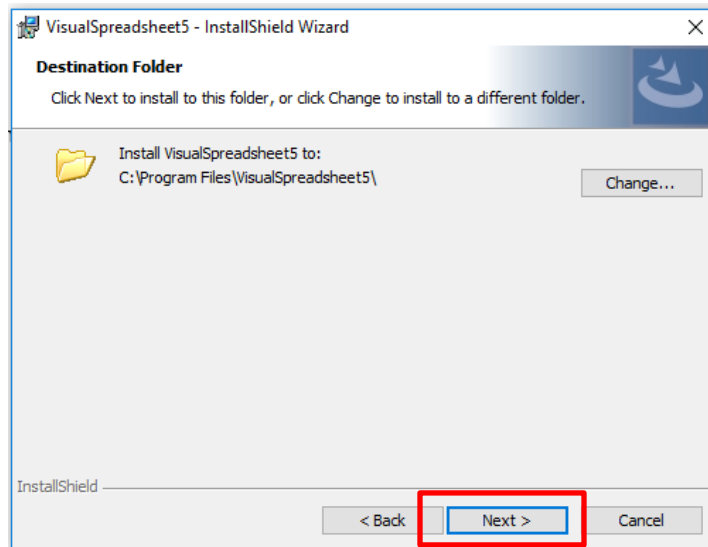


4. From there, you will see the Ready to Install the Program dialog, please **click INSTALL**.

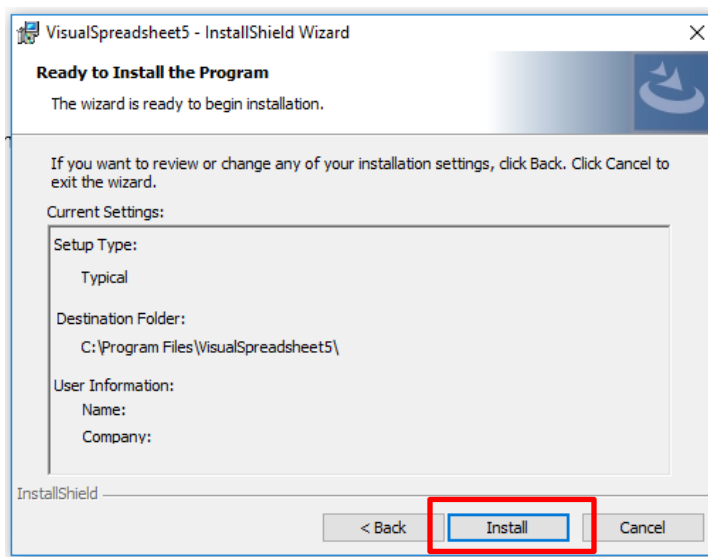


5. Please wait while VisualSpreadsheet 5 installs onto your system. Once it is complete, you will see the InstallShield Wizard Completed dialogue, please **click FINISH**.

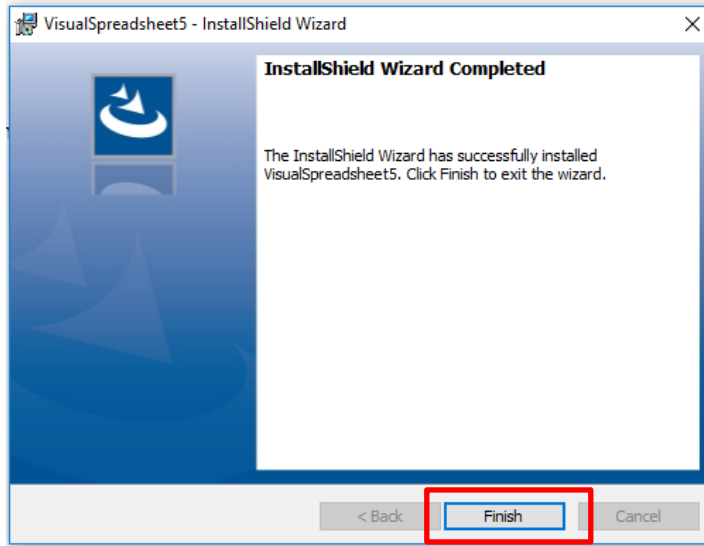


6. Please **RE-START** your system. This must be a full re-start of your system as this will initiate the installation of your new database.
7. Once your system starts back up, you will now be prompted to install the database which utilizes MariaDB software. When you see the Welcome to the MariaDB 10.2 (x64) Setup Wizard, please **click NEXT**.

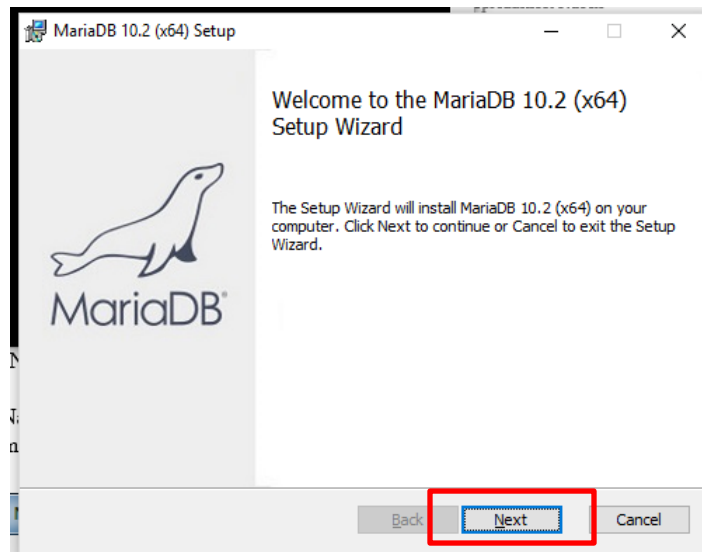

- You will then see the End-User License Agreement, once you have read the agreement, please **check the box** that says “I accept the terms in the License Agreement” and **click NEXT**.

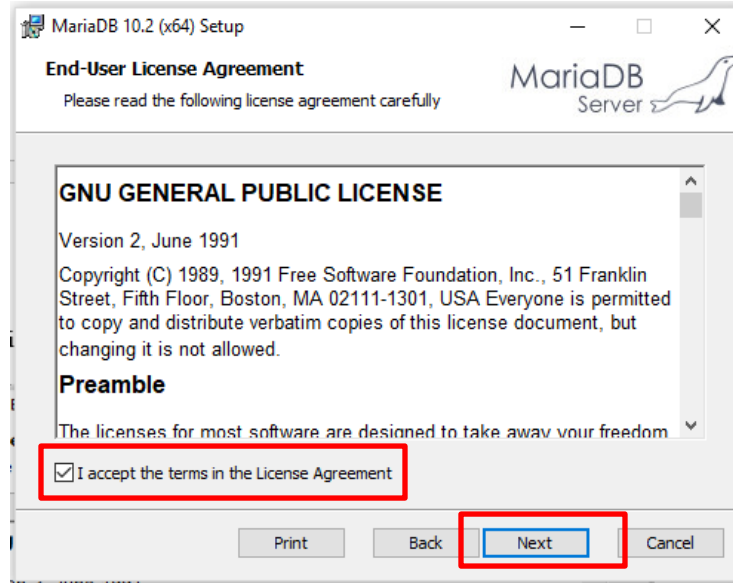


- The Customer Setup dialogue will then pop-up, **click NEXT**. *DO NOT* change any settings on this screen.

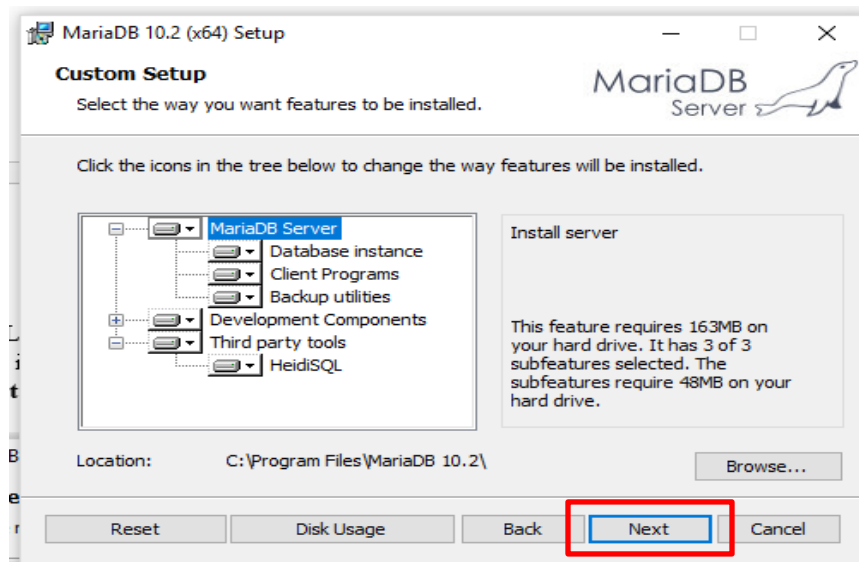


- You will then be prompted to set a root password in the Default instance properties dialogue. This root password must be set by the installer, we recommend using 8-12 characters.


Once you enter your root password, confirm and **click NEXT**. Please remember or write this password down, you will need it later to complete the installation! We recommend using the Serial number backwards followed by flowcam (for example SN10502: **20501flowcam**)

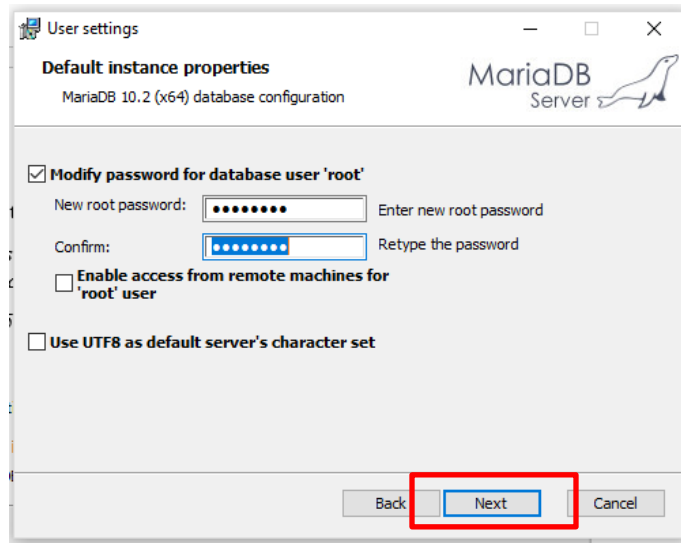


- Once your root password is confirmed, the Default instance properties dialogue will pop-up, please **click NEXT**. *DO NOT* change any properties or settings in this box.

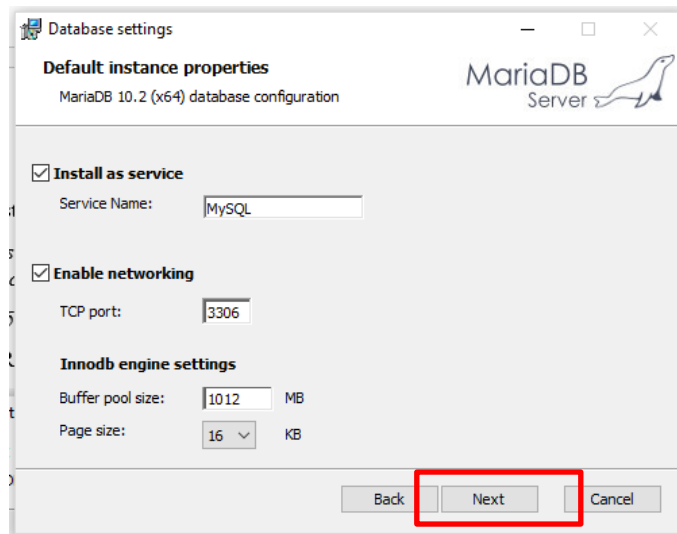


12. The MariaDB 10.2 Setup dialogue will now pop-up, please leave this box **unchecked** and **click NEXT**.

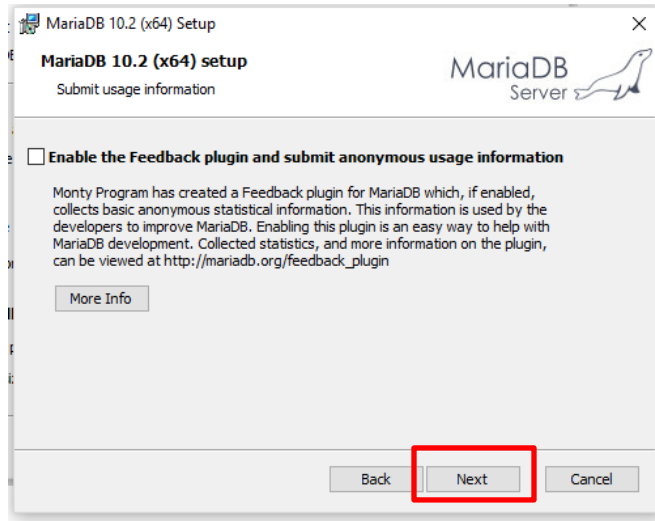


13. You will now see the Ready to install MariaDB 10.2 dialogue, please **click INSTALL**.

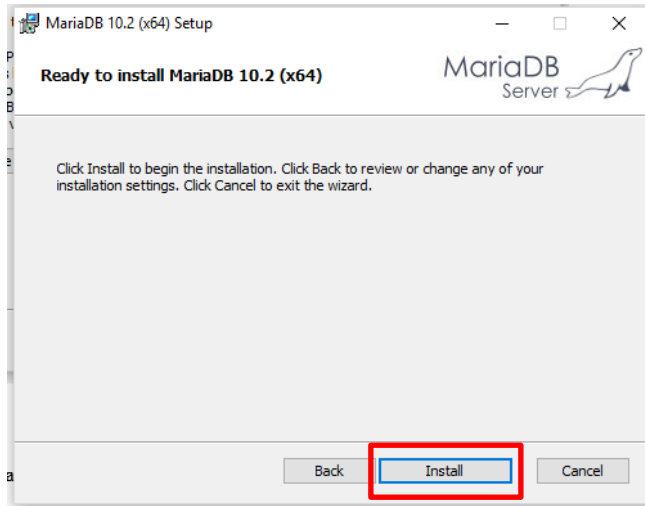

14. Installation will now begin and you will see a status bar, please *DO NOT* click anything in this box and allow the installation to complete.

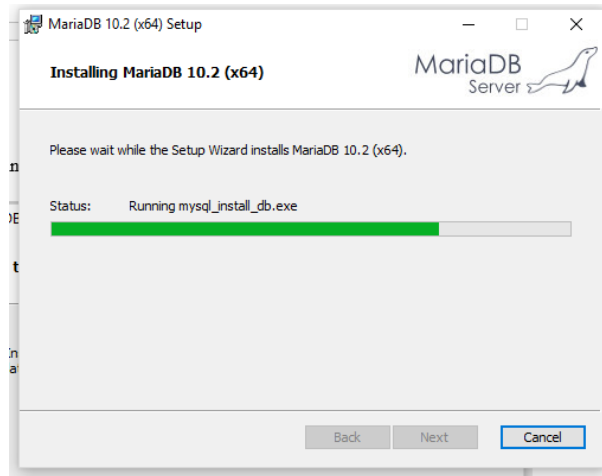

15. Once the installation is complete, you will see the Completed the MariaDB 10.2 Setup Wizard dialogue, please **click FINISH**. Your system will now do an automatic restart to finalize the installation of your new VisualSpreadsheet 5 database.

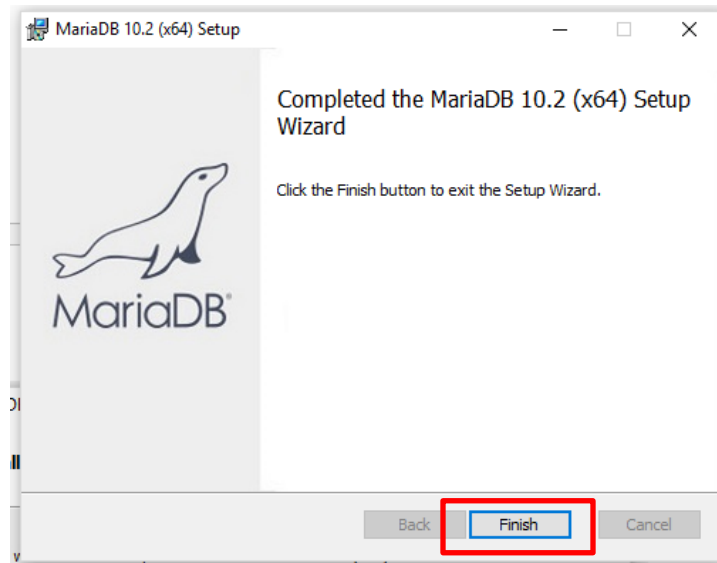


16. Open VisualSpreadsheet 5 by double clicking the icon on your desktop.





17. The Database Not Found dialogue box will now appear. **Please enter the following information:**

- a. Company: Your organization's name
- b. Password: This is the root password you created in Step 10
- c. Machine Name: If your FlowCam or PC has a unique name you can enter that here, alternatively you can use FlowCam followed by your serial number (i.e. FlowCam10502).
- d. Serial Number: This is the serial number of your FlowCam and can be found on the sticker on the right to where you install your flow cells.
- e. **Click OK**

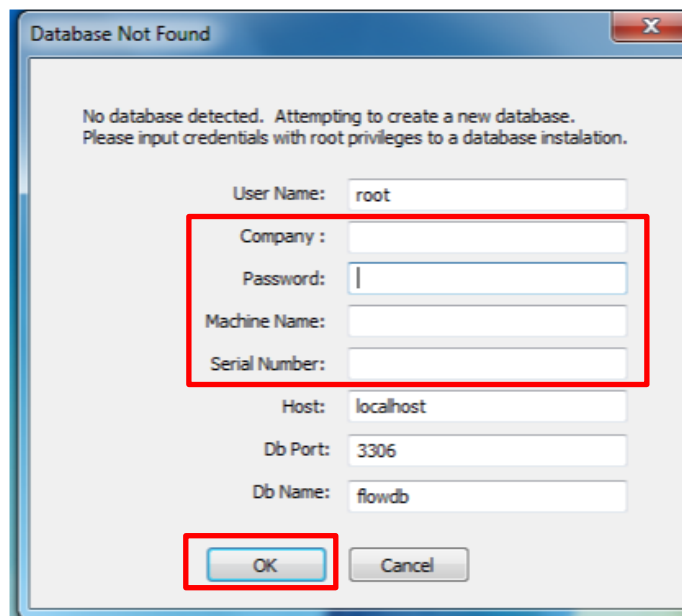

Database Not Found

No database detected. Attempting to create a new database.  
Please input credentials with root privileges to a database installation.

User Name: root

Company :

Password:

Machine Name:

Serial Number:

Host: localhost

Db Port: 3306

Db Name: flowdb

OK Cancel

VisualSpreadsheet will now open to the main page and the installation of VisualSpreadsheet 5 is complete.

\*Please note: If for some reason the installation does not proceed as expected, check in with your IT Administrator to assist you. Anti-virus software can prevent (eg Symantec) either MariaDB (database software) or the VisualSpreadsheet software installer to proceed properly. Your IT Administrator can temporarily disable any anti-virus software, so that VisualSpreadsheet can install properly.